

Scout[®] Pro Portable Electronic Balances



The OHAUS Scout® Pro – Hands Down The Industry's Best Portable Balance!

Designed for use in laboratory, industrial or education applications, the OHAUS Scout® Pro features easy-to-use two-button operation, a high-contrast LCD display, multiple weighing units, four application modes, and the option of either RS232 or USB connectivity, the Scout Pro is the portable balance for your needs!

- **Maximum Protection** Scout Pro is designed with mechanical and software protection against shock and overloading which can cause damage to the weighing cell. An integrated shipping lock protects the weighing cell during transportation or storage
- **RS232 or USB Connectivity** Scout Pro interface accessories are easy to install at anytime, even after purchase. The data cable is integrated in the kit, eliminating the need for additional cables
- Multiple Application Modes Scout Pro features Enhanced Parts Counting, Percent Weighing, Totalization and Display Hold

Scout Pro Portable Electronic Balancesunt

Built-In Application Modes:

Parts Counting — The Scout Pro is designed for counting parts for inventory control, parts bagging, receiving, dispensing and ticket counting

- Select from a sample size of 5, 10, 20 or 50 pieces while in counting mode with a quick button press
- The Scout Pro features user-enabled automatic recalculation of the average piece weight – instead of counting 100 parts for a sample, count 10 and let the Scout Pro count the rest of the sample! For applications where automatic recalculation is not desired, the easy-to-use menu system allows this feature to be disabled
- The Scout Pro can also be configured for situations where only the counting feature is required

Percent Weighing — Use the Percent Weighing mode to compare samples for compounding, formulation, or quick weight checking. Place your reference weight on the platform, enter the weight into memory, and the display will show the percentage of the sample weight on the platform compared to the reference weight.


Totalization— Use the Scout Pro's memory function to repeatedly sum the weight on the platform for compounding and formulation. Unlike other balances, the Scout Pro allows totalization up to the limit of the display!

Display Hold — The Display Hold mode allows you to retain the weight of an over-sized package on the Scout Pro's display. Great for postal and shipping applications!

User-Configurable — The Scout Pro allows you to select active weighing units and modes using the two button keypad. Scout Pro can be quickly configured using an on-display menu system – there are no cryptic codes to input! A built-in lockswitch prevents reconfiguration during use.

New Design for Portable or Benchtop Use



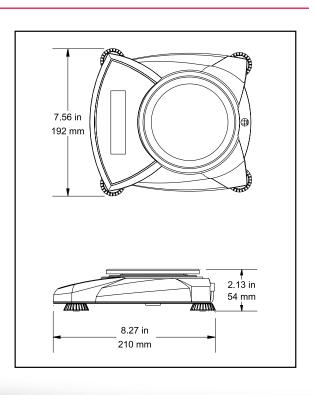

OHAUS Scout® Pro: Models and Weighing Units for Every Application

- · Parts Counting and Bagging
- Animal/Egg Weighing
- Feed Dispensing
- Medical Output Studies
- Crop and Seed Analysis
- Paint, Ink, Dye and Glaze Formulation
- Compound Preparation and Packaging
- Dental Adhesive and Compound Formulation
- Education Experiments
- Field Studies
- Parts Inventory, Quality Control and Receiving
- Filter Analysis
- Testing Kits
- Sample Gathering and Field Testing
- Food Portioning and Packaging
- Catering and Restaurant Applications
- Diet and Portion Control
- Jewelry Stores and Pawn Shops
- Injection Molding Sampling
- Shipping and Receiving
- Racquet and Golf Club Balancing
- Ticket Counting

The Scout Pro features the most common weighing units required for use in the lab, on the shop floor, or in the field:

Grams

Ounces

Troy Ounces

Kilograms

Pounds

Pennyweights

Combined Pound:Ounce display

Scout Pro Unit Availability

	60000	60.400	60400	60404	60404	600004	604664	604004	604000
Model	SP202	SP402	SP602	SP401	SP601	SP2001	SP4001	SP6001	SP6000
Capacity	200g	400g	600g	400g	600g	2000g	4000g	6000g	6000g
Readability	0.01g	0.01g	0.01g	0.1g	0.1g	0.1g	0.1g	0.1g	1g
g	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes
kg	-	-	-	-	-	Yes	Yes	Yes	Yes
OZ	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes
lb	-	-	Yes	_	Yes	Yes	Yes	Yes	Yes
lb:oz	-	ı	Yes	İ	Yes	Yes	Yes	Yes	Yes
ozt	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes
dwt	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes	Yes



using the included lockswitch

Scout[®] **Pro** Portable Electronic Balances

Model	SP202	SP402	SP602	SP401	SP601	SP2001	SP4001	SP6001	SP6000		
Capacity (g)	200	400	600	400	600	2000	4000	6000	6000		
Readability (g)		0.01			0.1						
Repeatability (Std. Dev.) (g)		0.01			1						
Linearity (g)	±0	.01	±0.02		±	0.1		±0.2	±1		
Capacity x Readability (kg)	N/A	N/A	N/A	N/A	N/A	2.0000 x 0.0001	4.0000 x 0.0001	6.0000 x 0.0001	6.000 x 0.001		
Capacity x Readability (oz)	7.0550 x 0.0005	14.1095 x 0.0005	21.645 x 0.0005	14.110 x 0.005	21.165 x 0.005	70.550 x 0.005	141.095 x 0.005	211.645 x 0.005	211.65 x 0.05		
Capacity x Readability (lb)	N/A	N/A	1.32280 x 0.00005	N/A	1.3230 x 0.0005	4.4090 x 0.0005	8.8185 x 0.0005	13.2280 x 0.0005	13.230 x 0.005		
Capacity x Readability (lb:oz)	N/A	N/A	1lb:5.165 oz x 0.001 oz	N/A	1lb:5.165 oz x 0.01 oz	4lb:6.550 oz x 0.01 oz	8lb:13.095 oz x 0.01 oz	13lb:3.65 oz x 0.01 oz	13lb:3.65 oz x 0.05 oz		
Capacity x Readability (ozt)	6.4300 x 0.0005	12.8605 x 0.0005	19.2905 x 0.0005	12.860 x 0.005	19.290 x 0.005	64.300 x 0.005	128.605 x 0.005	192.905 x 0.005	192.90 x 0.05		
Capacity x Readability (dwt)	128.60 x 0.01	257.21 x 0.01	385.81 x 0.01	257.2 x 0.1	385.8 x 0.1	1286 x 0.1	2572.1 x 0.1	3858.1 x 0.1	3858 x 1		
Application Modes	Parts Counting with Average Piece Weigh Recalculation, Percentage Weighing, Totalization, Display Hold										
Tare Range	To Capacity By Subtraction										
Over Range Capacity	Capacity +90d										
Stabilization Time	3 seconds										
Operating Temp. Range	50° to 104°F / 10° to 40°C										
Power Requirements	AC Adapter (included) or 4 AA Batteries (not included)										
Calibration Type	Digital Calibration from Keypad										
Calibration Mass (Full Span)	200 g	400 g	600 g	400 g	600 g	2000 g	4000 g	6000 g	6000 g		
Calibration Mass (Linear)	100 g	200 g	300 g	200 g	300 g	1000 g	2000 g	3000 g	3000 g		
Display	High			Contrast LCD with 6 digits 0.6in / 15mm high							
Pan Size (in/cm)	4.7 / 12 diameter		6.5 x 5.5 / 16.5 x 14.2								
Dimensions WxHxD (in/cm)				7.56 x 2.13 x 8.27 / 19.2 x 5.4 x 21.0							
Net Weight (lb/kg)	1.5 / 0.7			1.8 / 0.8							
Shipping Dimensions WxHxD	10 x 4 x 12 in / 25.4 x 10.2 x 30.5 cm										
Shipping Weight (lb/kg)		4 / 1.8									

Standard Features/Equipment

Removable stainless steel weighing platform, high contrast LCD display, lockswitch, integral security bracket, integral shipping lock, integral weigh below hook, sealed front panel and molded spill ring, stability indicator, overload and underload indicators, low battery indicator, auto shut-off, AC adapter, span calibration mass included on 200g, 400g and 600g models.

Accessories	OHAUS Number
Dot Matrix Printer	80252042
Adapter, RS232 Scout Pro Interface to the Dot Matrix Printer	80251934
Paper for Dot Matrix Printer	80251932
Ink Ribbon for Dot Matrix Printer	80251933
Thermal Printer	80251992
Paper for Thermal Printer	80251931
Adapter, RS232 Scout Pro Interface to the Thermal Printer	80251930
OHAUS Collect Software	80500746
TAL Software WinWedge	SW12W
RS232 Interface (cable Included) for SP, TA, NAV	
USB Interface (cable Included) for SP	71147377
Auxiliary Display Kit for Scout Pro	TAD7
Security Device for Scout Pro	76288-01
Scout Pro In-Use Cover, Set of 3	71168909
Hard Shell Carry Case for Scout Pro	77256-01

Industry Leading Quality and Support

All OHAUS Scout Pro balances are manufactured under an ISO 9001:2008 Registered Quality Management System. Our rugged construction and stringent quality control have been hallmarks of all Ohaus products for over a century.

OHAUS CORPORATION

* 7 Campus Drive Suite 310 Parsippany, NJ 07054 USA

Tel: 800.672.7722 973.377.9000 Fax: 973.944.7177

www.ohaus.com

With offices throughout Europe, Asia, and Latin America

*ISO 9001:2008

Registered Quality Management System



80774282_F © Copyright Ohaus Corporation